PWT1-PWT2 WASTE REFRIGERATOR USER MANUAL

English: PWT1-PWT2 Waste Refrigerator User Manual (Pages 1–17)

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1. Safety;

- Before making connections, check that the supply voltage is correct.
- The device should not be exposed to water or moisture. Protect the device from sudden temperature changes under high atmospheric humidity. **- Warning:** Before any maintenance, remember to disconnect all electrical connections.

1.1. General Information;

provides an overview of all important safety issues. Each section contains precise instructions for avoiding the hazards indicated by the above-mentioned warning signs. Following all important safety instructions ensures optimum protection against all hazards and ensures trouble-free operation. This device

has been designed using state-of-the-art technology . However, if the device is used incorrectly or for its intended purpose, hazards can occur

. To protect yourself from hazards, it is important to understand the contents of the user manual to avoid errors and operate the device safely and without errors. To avoid hazards and ensure optimum performance, no modifications to the device should be made that are not expressly approved by the manufacturer

1.2. Introduction: Please take a few minutes to read this manual carefully.

Proper maintenance and operation of this device will ensure the best possible performance from your product.

1.3. Package Contents;

1.3.1. Packaging, Storage, Disposal of Electrical and Electronic Products;

Before installation, completely remove the outer and inner packaging materials from the unit. Check that the device is complete and if any parts are missing, contact your customer representative immediately . When disposing of packaging, comply with the regulations in force in your country. Recycle reusable packaging materials . All packaging materials used are environmentally friendly. They can be stored without any risk or incinerated.

at the facility .

Correct Disposal of This Product (Waste Electrical and Electronic Equipment) (Applicable in the European Union and other European countries with separate collection systems)

This symbol shown on the product or its literature indicates that it should not be disposed of with other household waste at the end of its life.

To prevent potential environmental or human health harm from uncontrolled waste disposal, please separate it from other types of waste and recycle it responsibly to promote the sustainable reuse of material resources.

This product should not be mixed with other commercial waste for disposal and should be returned to a waste collection center.

1.4 Installation;

Note: If the device has not been stored or transported in an upright position, allow it to stand upright for approximately 12 hours before operating.

1- Remove the device from its packaging and ensure that all protective covering materials have been removed from all surfaces.

2- Place the device on a flat and solid surface.

1.5 DOCUMENTS;

The operating system of the product is produced and documented in accordance with European Standards (CE).

1.6 WARRANTY;

A new product is under warranty for 2 years.

Caution: Before performing any intervention on the product, the User and Technician must read this user manual carefully. Otherwise, the product may be voided from warranty.

1.7. Safety Instructions;

Information on occupational health and safety is based on European Union regulations valid at the time the device was manufactured.

If the device is used commercially, the user is responsible for ensuring that the occupational health and safety measures specified comply with the latest applicable regulations and that new specifications are adhered to throughout the entire life of the device.

Outside the European Union, occupational health and safety laws and local regulations applicable in the location where the device is installed must be observed. In addition to the occupational safety instructions in the user manual, general safety and accident prevention regulations relevant to the device's intended use, as well as environmental protection regulations, must be observed

⚡ Electrical Safety:

The device should only be used with a grounded socket.

It should not be used with an extension cord or multiple socket.

The device should not be plugged or unplugged with wet hands.

💧 Usage Safety:

The appliance should be used for brewing purposes only.

The ventilation openings of the device should not be blocked.

🔥 Fire and Explosion Risk:

The device should not be brought into contact with flammable and combustible materials.

Alcohol-based liquids should not be used for cleaning.

🔧 Care and Intervention:

User intervention is prohibited on the internal parts. In case of malfunction, contact an authorized service center.

The device should be turned off and unplugged before cleaning.

1. Product Description;

1.1 Information Usage ;

This manual is for the device It contains information on installation, operation, and maintenance . In addition to the information contained in this manual , the user must comply with all local health and safety controls and generally applicable safety regulations. The safe and correct use of the device covers the safety instructions and operating instructions contained in this manual. The user manual is part of the product and should be kept near the equipment and kept in a place easily accessible by personnel who will perform

Installation, maintenance, and cleaning. 2.1 General Information Description; This section

provides an overview of all important safety issues. Each section contains precise instructions for avoiding the hazards indicated by the use of the above-mentioned warning signs. Compliance with all important safety instructions ensures optimum protection against all hazards and ensures trouble-free operation. This device has been designed in accordance with the latest technology . However, if the device is used incorrectly and for its intended purpose, hazards can occur . To protect yourself from hazards, it is important to know the contents of the user manual to avoid mistakes and to operate the device safely and without errors. To prevent hazards and ensure optimum performance, no modifications should be made to the device that are not expressly approved by the manufacturer.

2.2 Safety Instructions;

Information regarding occupational health and safety is based on European Union regulations valid on the date of production of the device. In case of commercial use of the device, the user is responsible for ensuring that the specified occupational health and safety measures comply with the latest valid regulations and that new specifications are adhered to throughout the life of the device.

Internal depth: 43cm / 53cm (GN 1/1 Capacity) – 35-40 kg/m³

Environmentally Friendly HCFC-Free Polyurethane Insulation – Digital Control Panel – Automatic Defrost with Electrical Resistance – Special Water Evaporation System – Special Door Hinge System – Cleanable

and Removable Magnetic Door Seal – Outdoor Temperature: + 43°C & 65% Humidity

2.3 Technical Safety; -Before installation,

check the unit for visible external damage. Do not commission and/or operate a damaged unit .

-The appliance may only be connected to a properly installed single socket outlet that has protection and grounding. -Never unplug the power cord by pulling on the cord itself. Pull by holding the plug end. - Make

sure that the power cord does not touch heat sources or sharp edges. The power cord must not hang over the edge of the work surface .

Make sure that no one can step on or trip over the cord .

-The power cord must not be bent or tangled. -Do not use the appliance if it is damaged or not working properly. -Do not use accessories or spare parts not recommended by the manufacturer. This may cause damage to the user or personal injury and will void the warranty. -Do not move or tilt the appliance while it is operating.

- Never plug in the appliance with wet hands or damp hands. - Never operate the appliance on a wet or damp surface. - In case of fire, unplug the appliance or disconnect the appliance from the power supply before extinguishing. - Do not allow the appliance to come into contact with water while it is switched on. - Remove the appliance from its packaging and destroy all packaging materials. - Place the appliance on a flat and solid surface. - Never place the appliance on a wet, damp or flammable surface.

- Do not install the appliance near electric stoves or other heat sources. - Connect

the power cord of the appliance only if it complies with the regulation.

WARNING: If the mains cable of the appliance is damaged, contact the manufacturer or

distributor before operating it. - This appliance contains the refrigerant

propane (R290). Propane is a colorless flammable gas. It has a low global warming potential. It has no ozone depletion potential and replaces R12, R22, R134 and others. -Make sure that no part of the cooling device is damaged when moving or installing the device. -In addition to the operating noise of the compressor, flow sounds can also be heard in the entire cooling circuit. Unfortunately, these effects are unavoidable, but they have no effect on the performance of the device.

-The device should not be connected to the mains supply via extension cords or socket bars. These do not provide the necessary safety and may create a risk of overheating. -The manufacturer cannot be held responsible for damage caused by the missing or broken grounding conductor e.g., electric shock. -The electrical safety of the device is only guaranteed if it is connected to a protective grounding system installed in accordance with the regulations . It is important that this basic and technical safety requirement is met.

3.1 Delivery Check; Please

check the delivery for completeness and shipping damage upon receipt. In case of visible damage, please report.

Indicate the extent of the damage along with the carrier's delivery invoice.

3.2 Packaging; Before installation, completely remove the outer and inner packaging material from the unit.

Comply with the regulations applicable in your country when disposing of the packaging. Recycle reusable packaging materials.

Please check that the device is complete and if any parts are missing, contact your customer representative immediately.

3.3 Storage;

Keep the packaging closed until installation, in accordance with the installation and storage markings on the exterior.

The following notes should be observed when storing packages: - Do

not store in the open air. - Store in a dry and dust-free place. - Do not allow contact with corrosive substances and do not expose to direct sunlight. - In case of long-term storage (> 3 months) Regularly check the condition of the packaging and parts.

4.1 Safety Instructions;

-Do not use acidic substances and ensure that no water enters the device.

-To protect against electric shock, prevent the unit, cables or plug from coming into contact with water or other liquids. -The device is not suitable for direct cleaning with pressurized water . Therefore, do not use pressurized water to clean the equipment!

4.2 Cleaning;

-Clean the device thoroughly after each use.

-Before cleaning, disconnect the device from the power supply (unplug it) and wait until it reaches room temperature. -Open the lid and remove the waste bucket. -Wipe the device with a damp cloth to prevent liquid from entering the controllers . Clean the inside and outside of the unit, the door and the gasket with a damp cloth and a non-acidic cleaning agent.

- Thoroughly dry all cleaned surfaces - Always use a soft cloth; do not use abrasive cloths that could scratch the unit and equipment. - Dry the washed parts with a dry cloth. - Leave the door open for a while to allow the interior to dry completely. - If necessary, remove the door gasket and clean it.

4.3 Commissioning;

- Before operating the device, wipe the interior with a damp cloth and then dry it. - Plug the power plug into a grounded socket installed in accordance with regulations. The mains voltage must comply with the value data on the type label. - Turn on the main switch. The digital thermostat displays the current temperature in the cooling cabinet and the circulation fan operates continuously. The device should be operated until the factory temperature setting is reached . - Only the internally mounted ramp and garbage disposal can be removed and installed for easy replacement of waste disposal coolers . -Make sure that the air inlet and outlet are open. For the cooling device to operate properly, the above-mentioned warnings must be followed and the device must be clear of obstacles. - If the device will not be used for a long time, you should unplug it . Care should be taken to leave the covers and doors open to prevent odor formation in the uncooled interior.

NOTE: When you start the device; -If the device does not cool: Check the digital display of the device, are there any indicators ?

-Check that the power plug of the device is properly plugged into the socket? -Make sure that there is voltage in the mains socket? -Are the cover and door of the device opened frequently or should not be overfilled for cooling ? -Check that the device cover or door is closed properly ? -Check that the device door gasket is correctly seated?

4.4 Maintenance and Safety Instructions; -Installation, maintenance, and repair work

may be performed by specialists authorized by the manufacturer . The user is not responsible for any improper installation, maintenance, or repairs.

-Regularly check the power cord for damage. Do not use the device if the cord is damaged. If the power cord is damaged, have it replaced by an authorized service center.

-In case of damage or malfunction, please contact your dealer or customer representative. -Repair and maintenance of the device must only be performed by authorized personnel using original spare parts and accessories.

Note: Do not attempt to repair the device yourself .

Model: Models	Internal Temper ature (Temper ature °C)	240 Lt. Trash Can 240 Lt. Waste Box	Measur ements Dimensi ons Cm.	Packing Dimensi ons Packing Dim. Cm.	Weight Weight Kg.	Stainless Steel Type	Insulation Thickness Insulation Thickness	Cooling Gas Cooling Gas	Voltage Voltage
PWT-1	-2/+8	1 Name.	109x98x 128	118x105 x145	125	304	50 mm	R290	220V / 50hz

PWT1-PWT2 Trash Refrigerator User Manual: This detailed manual contains important information for the effective and safe use. If you encounter any problems, always seek assistance from a qualified technician.

Model: Models	Internal Temper ature (Temper ature °C)	240 Lt. Trash Can 240 Lt. Waste Box	Measur ements Dimensi ons Cm.	Packing Dimensi ons Packing Dim. Cm.	Weight Weight Kg.	Stainless Steel Type	Insulation Thickness Insulation Thickness	Cooling Gas Cooling Gas	Voltage Voltage
PWT-2	-2/+8	2 Names.	174x98x 128	183x105 x145	175	304	50 mm	R290	220V / 50hz

PWT1-PWT2 Trash Refrigerator User Manual: This detailed manual contains important information for the effective and safe use. If you encounter any problems, always seek assistance from a qualified technician.

Using

digital ;

SET	the set value ; confirms the values in programming mode.
	(DEF) Starts Manual Defrost .
	(UP) : Shows the maximum measured temperature; in programming mode, it is used to change parameters and increase values.
	(DOWN) : Displays the minimum measured temperature; In programming mode, it is used to change parameters and decrease values.
	Used to turn off the device (onF = Off whereas) It is used.

KEY COMBINATIONS;

+	Locks and unlocks the keypad.
SET +	To enter programming mode
SET +	To exit programming mode

Use of LEDs

The function of each LED is described in the table below .

LED	MODE	FUNCTION
	OPEN	The compressor works.
	Flashing	Flashes Between Compressor Operations.
	OPEN	Defrost Works.
	Flashing	Drainage Time
	OPEN	Fans On
	Flashing	Fan Delay After Defrost
	OPEN	Alarm Status Available.
	OPEN	There is a continuous cycle.
	OPEN	Energy saving is achieved.

°C/°F	OPEN	Unit of Measurement
°C/°F	Flashing	Programming Phase

Seeing the Minimum Temperature

1. Press and release the ▼ key.
2. The message "Lo " is displayed , followed by the recorded minimum temperature.
3. Pressing the ▼ button again or waiting for 5 seconds returns to normal display.

Seeing Maximum Temperature

1. **HE** Press and release the key.
2. The " Hi " message is displayed, followed by the maximum recorded temperature.
3. Press the **O** key again or wait 5 seconds to return to normal display.

Resetting Recorded Maximum and Temperature Values

1. While the maximum or minimum temperature is displayed, press and hold the **SET** BUTTON for more than **3 seconds** (the "rSt" message will appear .)
2. The "rSt" message will flash and return to normal temperature display.

Displaying the Setting Value

1. Press the **SET** button : the setting value is displayed. 2.

Press the **SET** button or wait **5 seconds** to exit .

Changing the Setting Value

1. Press and hold the **SET** button for **2 seconds**.
2. The set value is displayed and the display unit flashes "C" or "F" .
3. ▼ or Adjust the setting value with the **O** key.
4. After setting the value, confirm the value by pressing the **SET** key.

Starting Manual Defrost

Defrosting starts when you press and hold the **DEF** button for **2 seconds** .

Locking the keys

1. ▲ Press and hold the + ▼ keys for **3 seconds**.
2. When the keypad is locked, the "**POF**" message will be displayed.
3. When a key is pressed for 3 seconds, the "**POF**" message will be displayed again.

Unlocking the keypad

1. ▲ Press and hold the + ▼ keys for 3 seconds. The "PON" message will appear on the screen and the key lock will be unlocked.

Continuous loop

defrosting continues , it is activated by pressing the ▲ key and holding it for **3 seconds**. The compressor is activated to maintain **the CSS** setpoint for the time set in the **CCT** parameter. **The cycle can be terminated before the end of the setpoint by pressing the same** activation key, the ▲ key, for 3 seconds.

MESSAGE	WHY	EXITS
"P1"	Room sensor error	Compressor output depends on the " Con " and "COF' " parameters .
"P2"	Evaporator probe failure	Defrost finish is timed.
"P3"	Third sensor error	There is no change in outputs
"P4"	Fourth sensor error	There is no change in outputs
"HA"	Maximum temperature alarm	There is no change in outputs
"LA"	Minimum temperature alarm	There is no change in outputs
"HA2"	Condenser high temperature	It depends on the "Ac2" parameter.
"LA2"	Condenser low temperature	" Bll " parameter.
"In"	Open door	Compressor according to parameter " rrd "
"EA"	External alarm	There is no change in outputs
"CA"	Serious external alarm (i1F=Bal)	All exits are closed
"CA"	Pressure automatic alarm (i1f=PAL)	All exits are closed

Alarm Correction

Sensor errors "P1," "P2," "P3 ," and "P4" become active a few seconds after an error is detected in the relevant sensor; the alarm signal is interrupted after a few seconds of normal sensor operation. It is beneficial to check the connections before replacing the sensor. Temperature alarms

"HA," "LA," "HA2 ," and "LA2" are interrupted immediately after the temperature returns to normal values.

" EA" and "CA " alarms (i1F=Bal) are interrupted after the digital input connection is disconnected.

" CA" Alarm (i1F=PAL) is interrupted by rebooting the device.

Other Messages

Pon	The keypad is unlocked.
PoF	The keypad is locked.
noP	programming mode: There is no parameter in Pr 1 step.
noA	On the display or at Dp 2,dp3,dp 4: The selected sensor is not in use.

EVCO Controller Usage and Installation Instructions;

1-General Warning :

Digital thermometer with relating to 30 per second One process not done time , digital thermometer self locks . First when pressed to the screen « Loc » article you will see . This open for any One press the button for 2 seconds printed hold . « UnL » text will be seen And from now on digital process to do possible will be .

Key team's for use opening (« Loc » text on the screen)

-Press any button for more than 2 seconds. The text " UnL " will appear on the screen. You can use the keys. For security reasons, if you don't press any keys for 30 seconds, the device will automatically enter key lock.

To perform any action, first remove the key lock

Viewing and changing the set value

- **SET** Press the button once quickly and then release your hand. The set value will appear on the screen and the LED will flash.
- or keys to set it to the required value within 20 seconds
- **SET** Press the button again or wait 20 seconds without pressing any buttons.

Note: The set value can be changed between the values assigned by the r1 and r2 parameters. Please refer to the manual for detailed information.

Defrost

defrosts every 8 hours according to factory settings .

Note: This time period starts as soon as the device is powered on.

Starting Manual Defrost

- Press the DEFROST button and hold it for 3 seconds.
- The defrost cycle will start immediately and the LED will light. According to the factory setting, this defrost will last 20 minutes.

Note: A manual defrost cycle always resets the defrost counter; a new defrost begins after 8 hours (or after the assigned d0 time).

If defrost values are changed while the device is being initially configured, performing a manual defrost once after the initial setup resets the current defrost cycle and allows defrosting to continue according to the newly assigned values .

TURNING THE DEVICE ON AND OFF

- You can switch the device to and from standby mode by pressing and holding the button for 4 seconds. The LED will flash while you hold the button, indicating that the operation is in progress. If no text appears on the screen and only the red LED in the lower right corner is lit, the device is off. Please turn the device on before any further operations.

IMPORTANT WARNINGS

Control panels provide users with various functions to prevent the spoilage of products stored in the refrigerator and to easily identify potential issues.

IMPORTANT WARNING: Turning off your product via the control panel only puts it into standby mode. To cut off power to the control panel, turn off the power switch on your product's electrical box or unplug it from the outlet.

-ALARMS

AL: Low temperature alarm. An icon will appear on the control panel indicating which probe is faulty, such as PR1, PR2 or PR3. Contact technical service.

AH: High temperature alarm. When loading your device, be sure to follow the instructions in the user manual. After ensuring the door is closed, your device should reach the desired temperature within 60 minutes while empty. Please contact authorized service.

PR1 : Cabin sensor error. Please contact authorized service.

PR2 : Evaporator sensor error. Please contact authorized service.

ID : Door switch alarm. If the door remains open for more than 1 minute, the alarm will sound. The alarm will go off when you close the door. If the alarm persists, check the door switch connections. If the alarm persists, contact authorized service.

IA : Power outage alarm. If the power supply to the system is cut off, the alarm will start. Contact authorized service.

COH : Indicates that the temperature detected by the condenser sensor has reached the first critical point. Please unplug the device and clean the condenser . Wait 30 minutes before restarting your device. If the alarm persists, please contact authorized service.

CSd : Indicates that the temperature detected by the condenser sensor has reached the upper limit. In this case, the compressor will be deactivated and the device's power supply will be cut off. Please check and clean the condenser . Wait 60 minutes before restarting the device. If the "COH" or " CSd " alarm occurs again, contact your authorized service center.

WARNING: If the intervention specified as a result of COH and CSd alarms is not performed, compressor malfunctions that may occur will be excluded from the warranty.

dFd : Indicates that the device has been defrosting for the maximum time , but frosting continues because it has not reached the desired defrost temperature. Please manually defrost the device again ; to do this, press and hold the defrost button on the display for 4 seconds. If the alarm persists, contact authorized service.

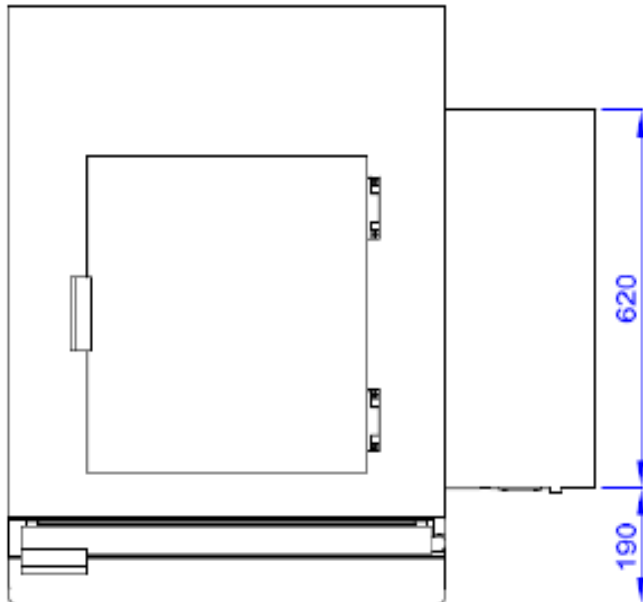
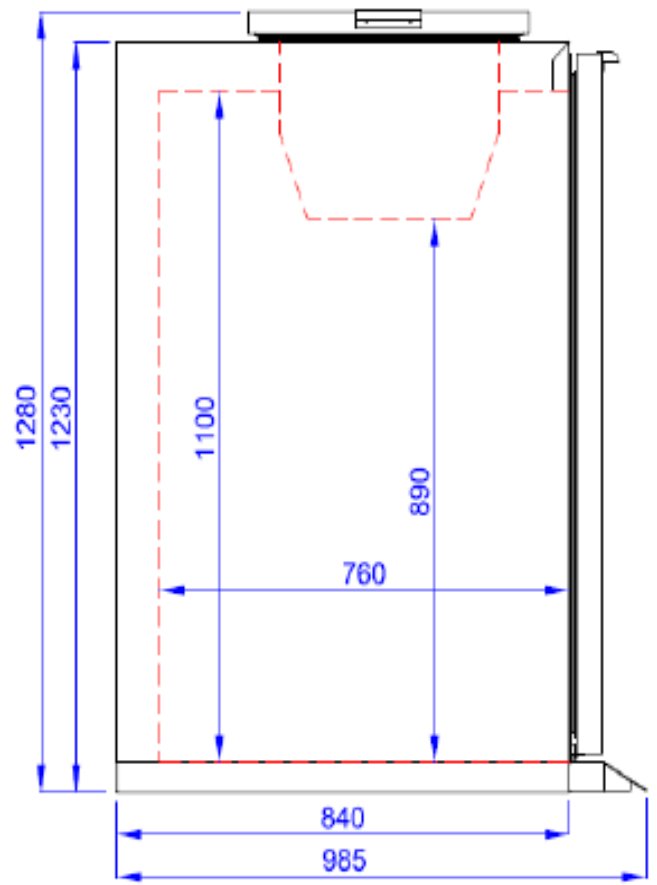
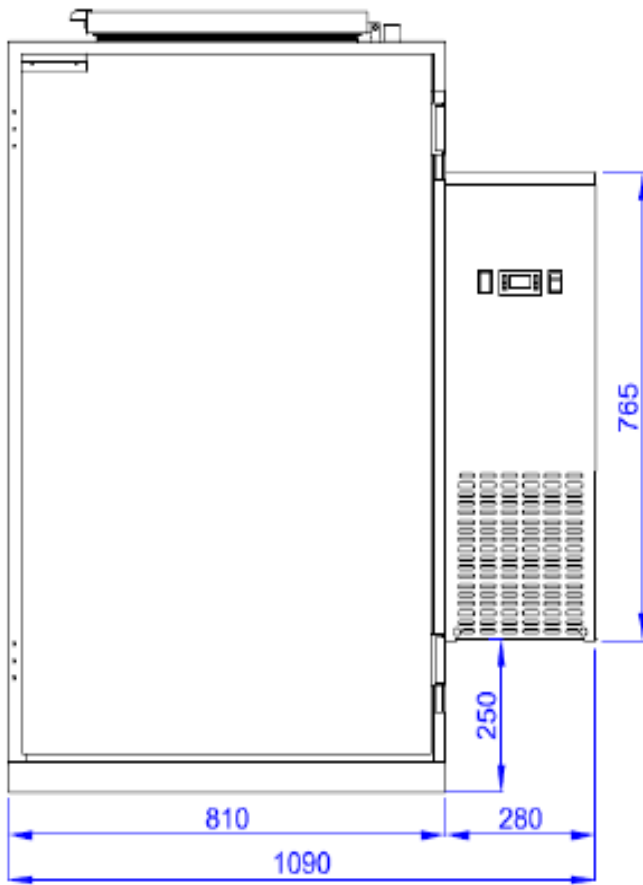
WARNING: DIGITAL CONTROL PANELS ARE SPECIFICALLY ADJUSTED IN OUR FACTORY TO ENSURE THE MOST CORRECT OPERATION OF THE PRODUCT. NEVER CHANGE ANY CONTROLLER PARAMETERS OTHER THAN INTERNAL TEMPERATURE ADJUSTMENT. OTHERWISE, SERIOUS DECREASES OR PROBLEMS MAY OCCUR IN THE DEVICE'S PERFORMANCE.

ALARM CODE	MEANING	WHY	WHAT NEEDS TO BE DONE
P1	Media probe reading failure	The ambient temperature sensor may be damaged or broken.	Call technical service.
P2	Evap probe reading failure	The evaporator temperature sensor may be damaged or broken.	Call technical service.
in	Door open alarm	The door does not close or the door switch is faulty.	Is there anything preventing the door from closing? If not, call technical service.
HA	High temperature alarm	an icing or cooling problem.	Manual defrost if there is frost You can enter defrost mode by pressing and holding the snowflake icon on the digital display for 3 seconds . If the problem persists, call technical service.

Changing the Cabinet Operating Temperature

DIXELL

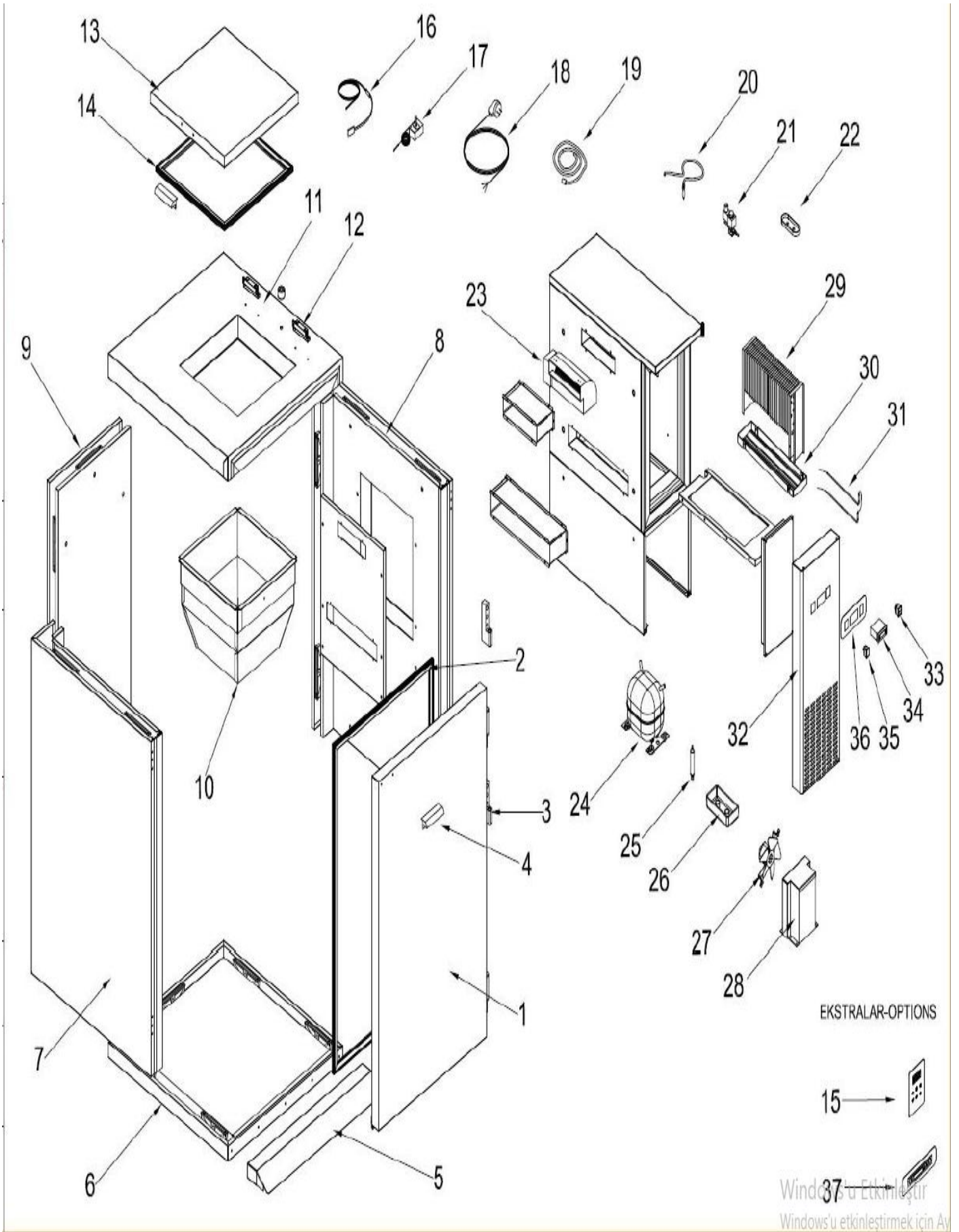
Press and hold the "SET" button for 3 seconds. When the "C" symbol in the upper right corner flashes, select the desired temperature with the "up and down arrow" buttons. Then press the "SET" button.

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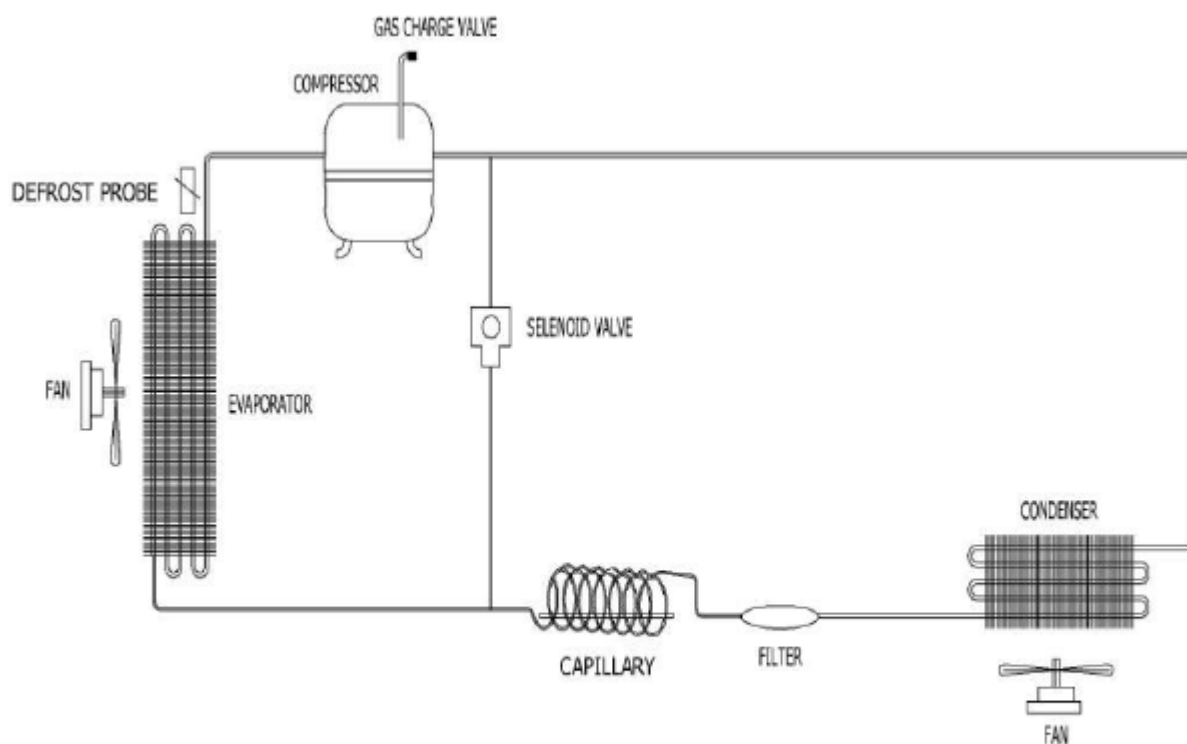
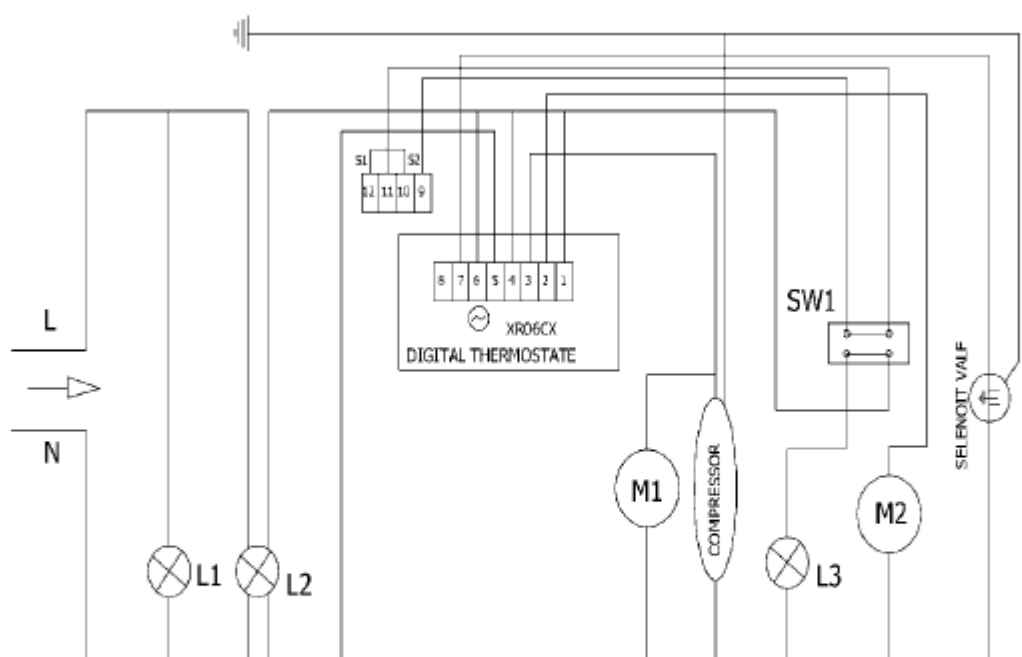
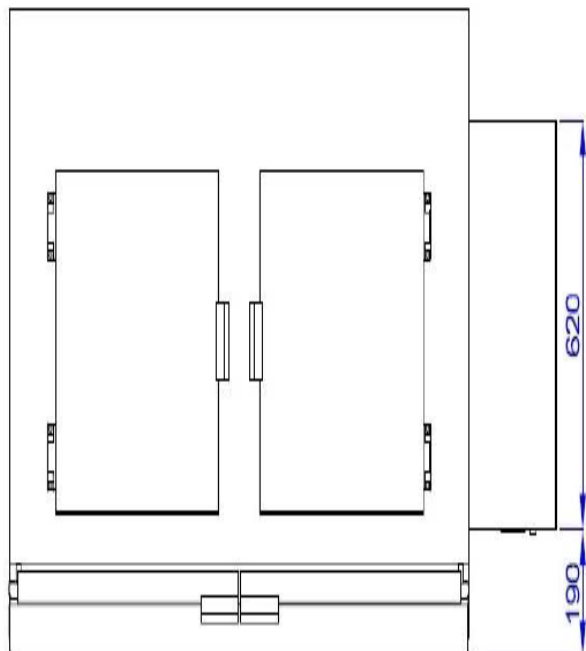
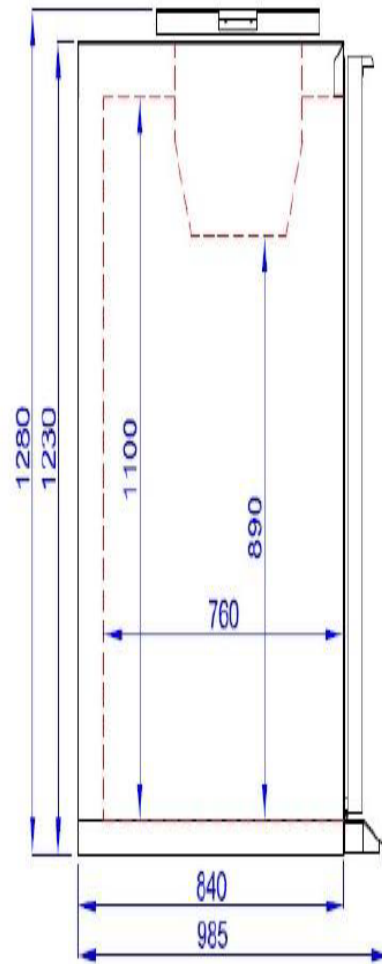
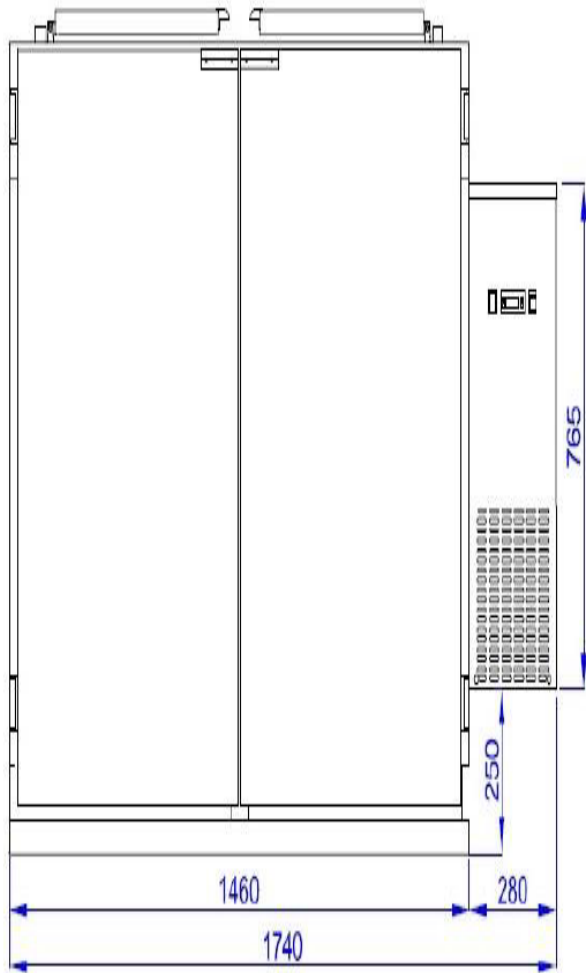
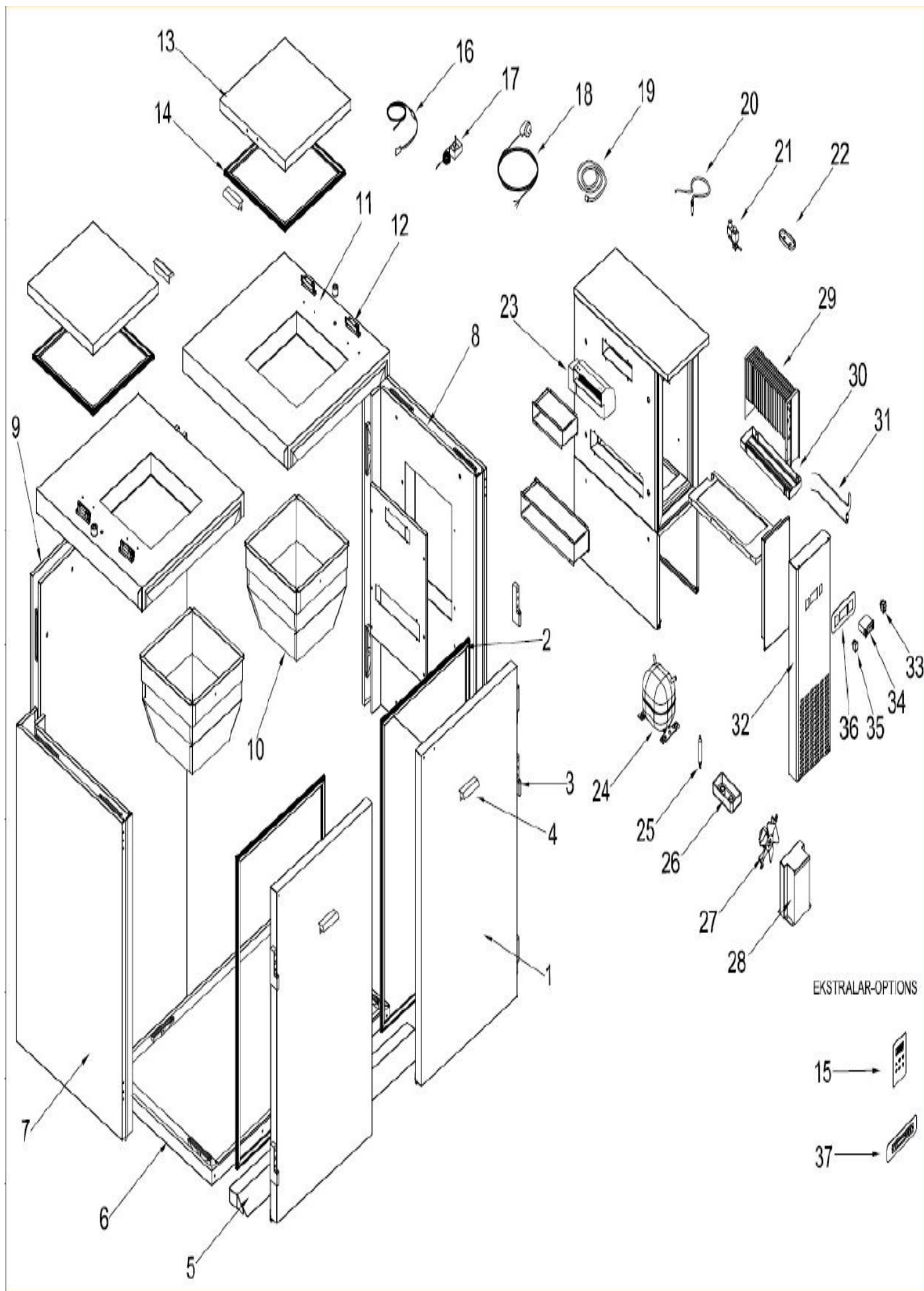
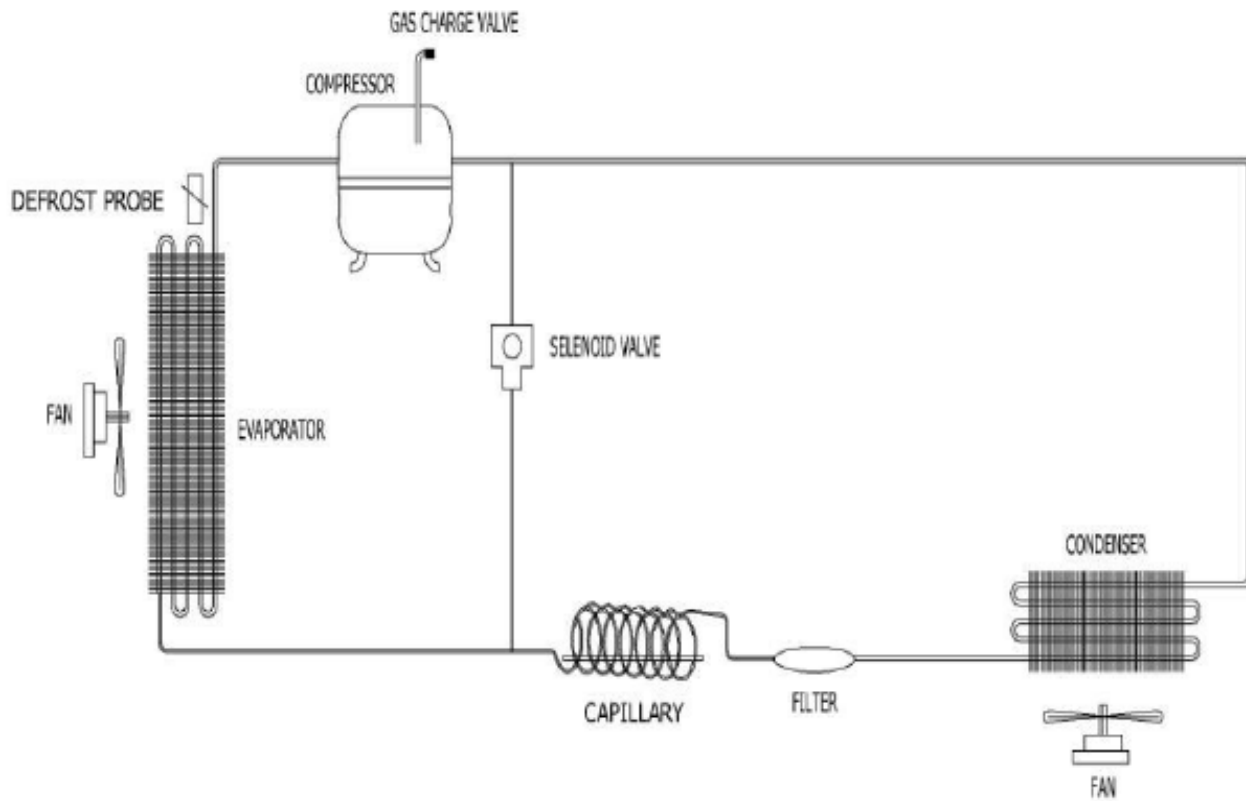
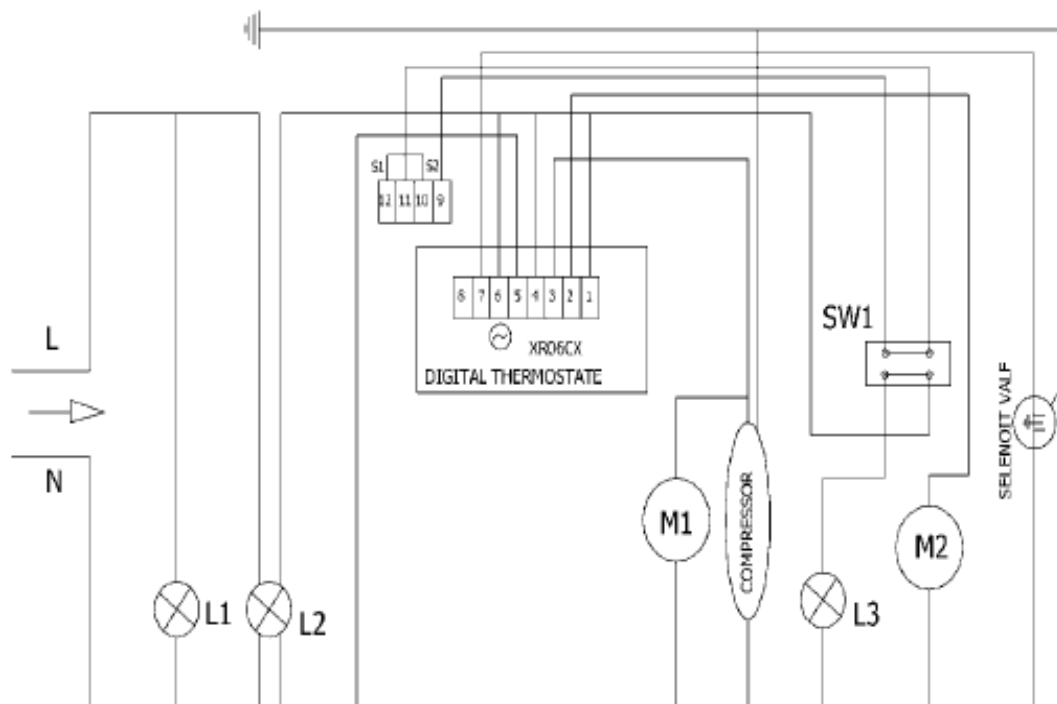
L1 : POWER LAMBASI / POWER LAMP
L2 : ACMA KAPAMA LAMBASI / ON OFF LIGHT
L3 : KANOPI LAMBASI / DEFROST LIGHT
SW1 :ON/OFF SWITCH

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